

# GUIDE | TAGARNO RULER APPLICATION

VERSION: 1.2 | 2018-02-01

Use the control box to access the setup menu, go to Ruler under the Applications menu point to turn the Ruler application on horizontally, vertically or both directions and get a ruler displayed on top of your live image on screen.

Turn the ruler off by once again accessing Ruler option under the Applications menu point and choose Off.



## System units

Choose to display metrics or imperial units when using the Ruler. Go to the System units option under the General settings menu point in the setup menu.



## Ruler color

Set the ruler color to your preferred color choice to separate it from your displayed object. Go to the Ruler color option under the General settings menu point in the setup menu.



Choose between 6 different colors: Yellow, blue, red, black, white and green.

## Manual focus setting

When the Ruler is chosen the system will automatically start in manual focus setting to focus on the part of the object furthest away.

If needed, manually adjust the focus settings and the height of the microscope to get the image in focus.

Please use "Optimal Working Distance" in Camera settings menu, before using Ruler



## Save image with or without graphics

It is also possible to save an image with the Ruler graphics by long-pressing the image button on the control box. Short-pressing the button will save the image without graphics.