


# GUIDE | TAGARNO WATERMARK APP

VERSION: 1.2 | 2018-02-01

Create an info field on captured images with an image prefix, time stamp, horizontal field of view, microscope model plus serial number and/or import and add a custom defined watermark.

When activated, the watermark will now appear on all captured images.

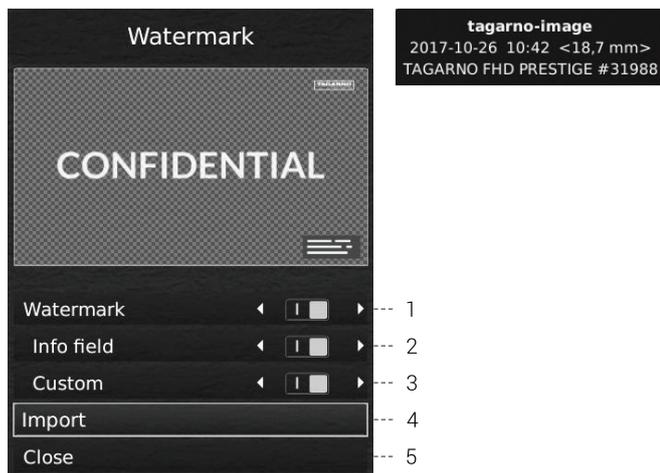

## 1. Enable or disable all settings

Choose to enable or disable all watermark settings at the same time.

## 2. Enable or disable info field

Enable or disable an info field, automatically stating an image prefix (default or as stated under Camera settings), time stamp, horizontal field of view plus microscope model and serial number.

## 3. Enable or disable a custom watermark

Choose to enable or disable the latest imported custom watermark.

## 4. Import a custom watermark

At this stage, you are able to import a custom watermark by inserting a USB memory stick into the microscope containing a watermark file.

The file must be a PNG type, measure 1920x1080px and be named watermark.png.

It is possible to import only one custom watermark at a time. This means that if you wish to change the current watermark, you need to follow the steps above to import a new watermark.

The application provides a range of options to control the watermark:

1. Enable or disable all settings
2. Enable or disable info field
3. Enable or disable a custom watermark
4. Import a custom watermark